

## **AET RTA Operating Instructions**

1. Open door by double clicking "Open Door" in the menu.
2. Lift door lever and slowly pull door open.
3. Load sample.
4. Close door and press door lever door down.
5. Close door by double clicking "Close Door" in the menu.
6. Select "Start Recipe" from menu.
7. Select the "Advanced" tab.
8. Select "Browse".
9. Select company or group folder in E:\Recipes.
10. Select recipe.
11. Select Execute.
12. After recipe is complete, select "Open Door" from the menu.
13. Remove sample.
14. Select "Close Door" from menu.
15. Fill out logbook.
16. Complete.